3. Beleuchtung

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EN

1. Viewing location

Before use, select a suitable location to set up the microscope. It is important to have sufficient lighting available (window, desk lamp).

2. Normal viewing

Place the object you wish to view (e.g. a stone) in the middle of the microscope table. The larger the object, the easier it is to fix it under the object clips (6). Look through the binocular lens (1) and adjust the distance of the eyepieces so as to obtain a clear, round image. You can sharpen the image by slowly turning the focusing wheel (7).

3. Light

Remove the two eyepieces (3) from the lens tubes (1) and turn the microscope around. The battery compartment is situated under the microscope's base.

(9). Press on the clasp (10) and open the battery compartment by lifting the cover (11) at an angle (while continuing to press on the clasp).

Place the batteries in the compartment so that the flat negative poles (-) press firmly against the spring terminals and the positive poles (+) touch the flat contacts. Close the battery compartment's cover (11) and turn the microscope back up.

Reinsert the eyepieces. You can now switch on and off the light (4) using the On/Off switch (8) at the base of the microscope.

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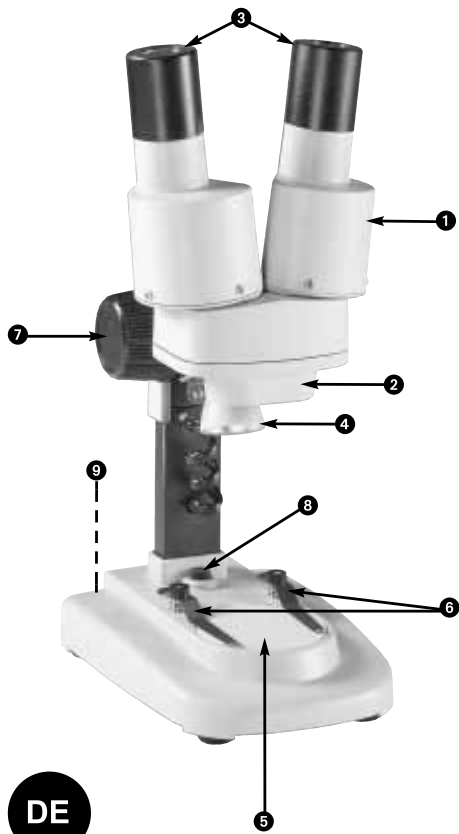
EN

Parts (Fig. 1)

- 1. Binocular view (with eyepieces)
- 2. Binocular lens
- 3. Eyepieces
- 4. Lighting (LED)
- 5. Microscope table (with plate for objects)
- 6. Clips for objects
- 7. Focus wheel
- 8. On/Off switch
- 9. Battery compartment
- 10. Battery compartment clasp
- 11. Battery compartment cover
- 12. Battery contacts (flat and spring terminals)

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