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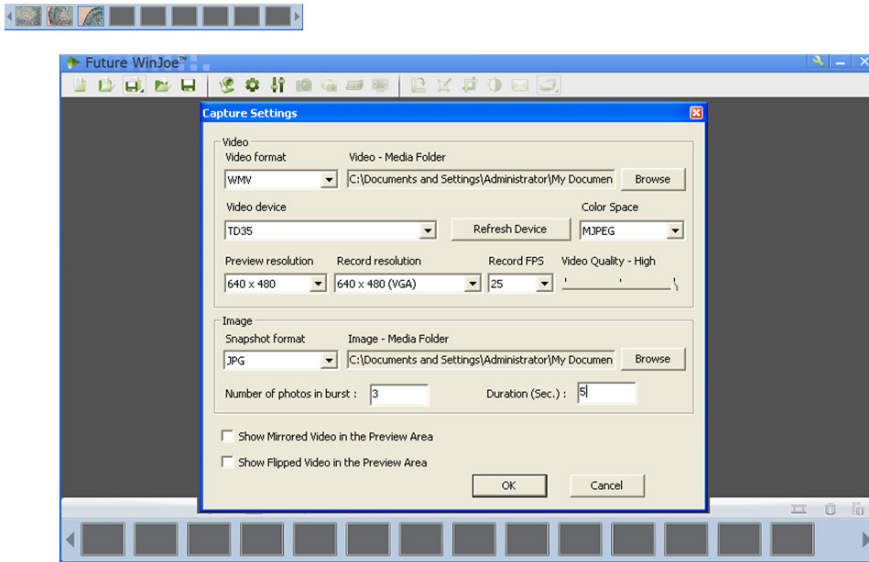
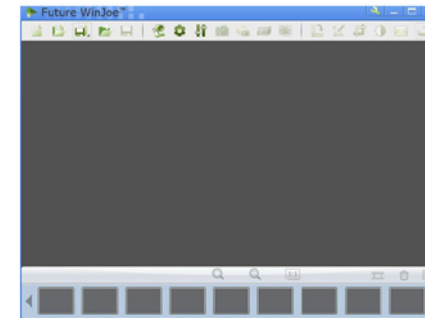
9. Click  to start video capturing, and click again to stop.

10. In order to view the pictures you have captured, you should click the pictures and videos in the list area below. After viewing, to click  to start another capture

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
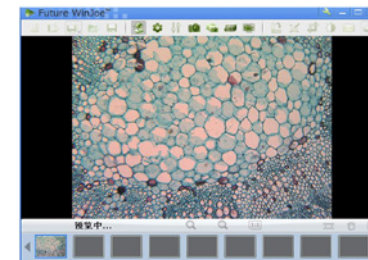
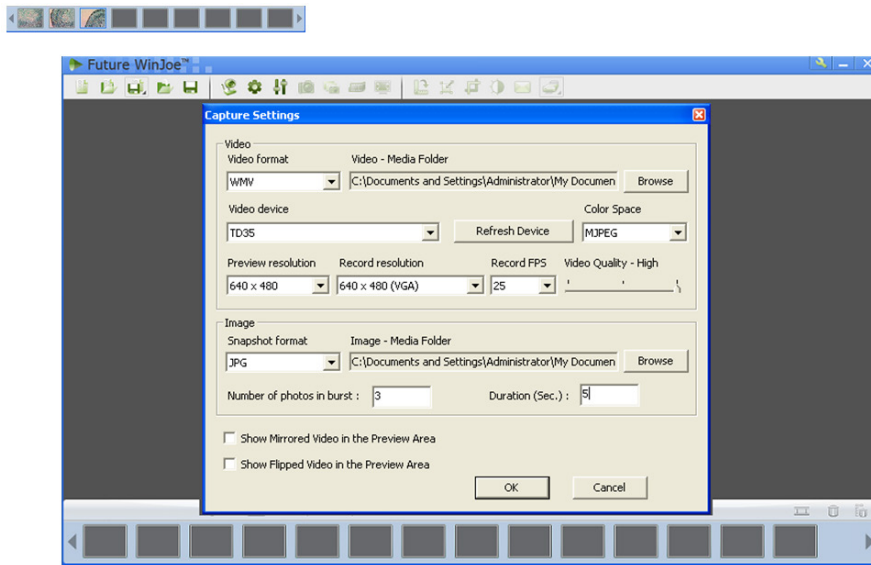
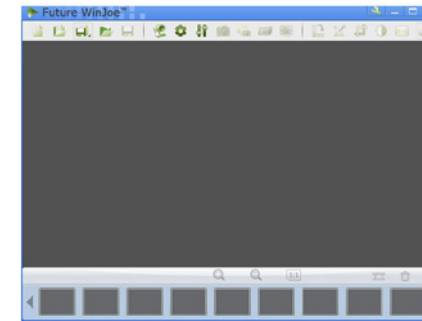
OPERATION OF FUTURE WIN JOE

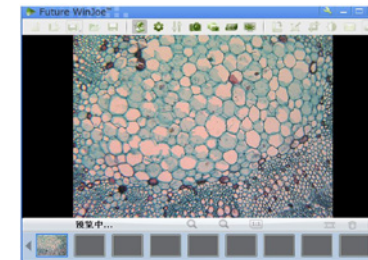1. Connect the digital device to the USB port,
2. Double click icon on desktop «  ».
3. There are two main parts in the interface: up-part for alive image showing and the down-part for list of images captured


4. Click»  « to let digital device capture images
5. In this manual, we display biologic microscope and digital eyepiece, please adjust the focus-nob of a microscope to get a clear image.


6. Before capture, please set the destination folder for saving image files.
7. Click»  «, to set the attributes of images shown, including brightness, contrast and so on.
8. By click»  «, you can get a table of “capture setting” and by changing the related value, and click  again to turn on capturing, you can view images with new specifications

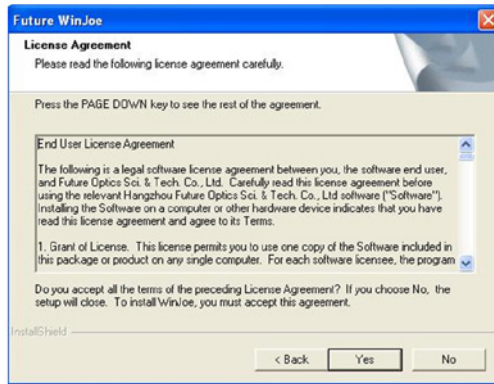
5. You must accept the user license agreement, so please click «yes» to continue.

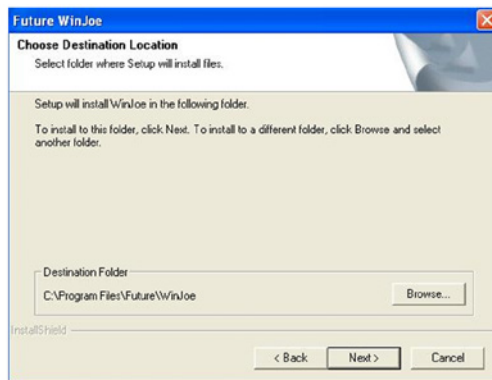6. Win Joe, the default location is «C:\Program Files\Future\WinJoe», of course, you can change it to the destination folder you want to click "Browse"

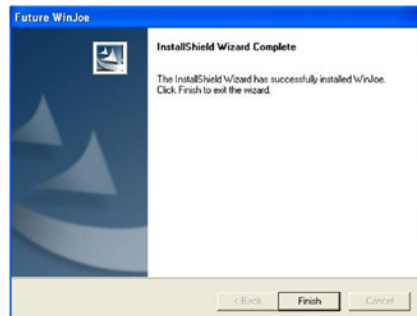7. Install shield wizard has been installed on your computer, click «finish» to exit.

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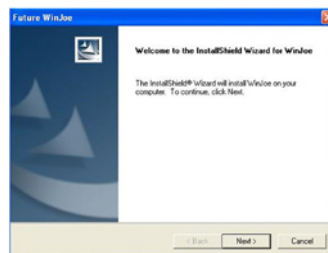
SOFTWARE FUTURE WIN JOE INSTALLIEREN

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DRIVER INSTALLATION

Please click «Install MD35 driver» to install the MD35's driver. Note: Please do not Insert the electronic eyepiece into the computer.

To connect Digital Eyepiece

Before connect the Digital Eyepiece to your PC, please confirm that the USB1.1 or USB2.0 interface on your PC is all right.

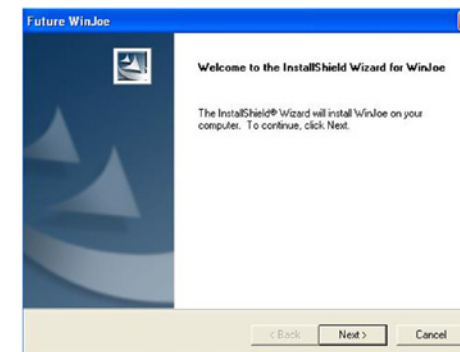
- Take out one optic eyepiece of the using Microscope, put in the digital eyepiece.
- Connect Digital Eyepiece to your PC
- “ Find a new hardware—” will appear on the computer screen. Wait for a moment, the electronic eyepiece can be work.

INSTALL SOFTWARE OF FUTURE WIN JOE

1. Put the CD into the CD-ROM, the files of CD will be shown on the screen automatically.
Notice : If the CD-ROM can not run automatically, please double click “My computer” then click the CD with right key or double-click “setup” under the folder of “software” of CD.
2. Click “Install Future Win Joe” option.
3. Select your language, then click “OK” to confirm

4. Come to the welcome wizard interface, click “next” to continue.

MD35 Digital Eyepiece for Microscope Installation Instructions

PLEASE CAREFULLY READ THIS MANUAL AND KEEP IT IN A SAFE PLACE.

INTRODUCTION

Congratulate for you owning the MD35 Digital Eyepiece. This Eyepiece has 0.35M pixel, it is suitable for any standard biological microscope, stereo microscope or metallography microscope. The Eyepiece can change images into digital signal, and send it to your PC, so that you can obtain, analyze the digital image and communicate with other people. Also the Eyepiece can work with other tools, such as projection machine, CD rewriter, and notebook PC, and it will make your work easy, interesting, and effective.

PACKAGE CONTENTS

When first opening the packing box, note carefully the following parts:
·MD35 Digital Eyepiece with usb cable · CD-ROM · Adaptor ring ·

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